
**Operating instructions**  
**Bedienungsanleitung**  
**Mode d'emploi**

**METTLER TOLEDO**

**Micro draft shield for AT21/AX26**  
**Mikro-Windschutz zu AT21/AX26**  
**Micro pare-brise pour balances AT21/AX26**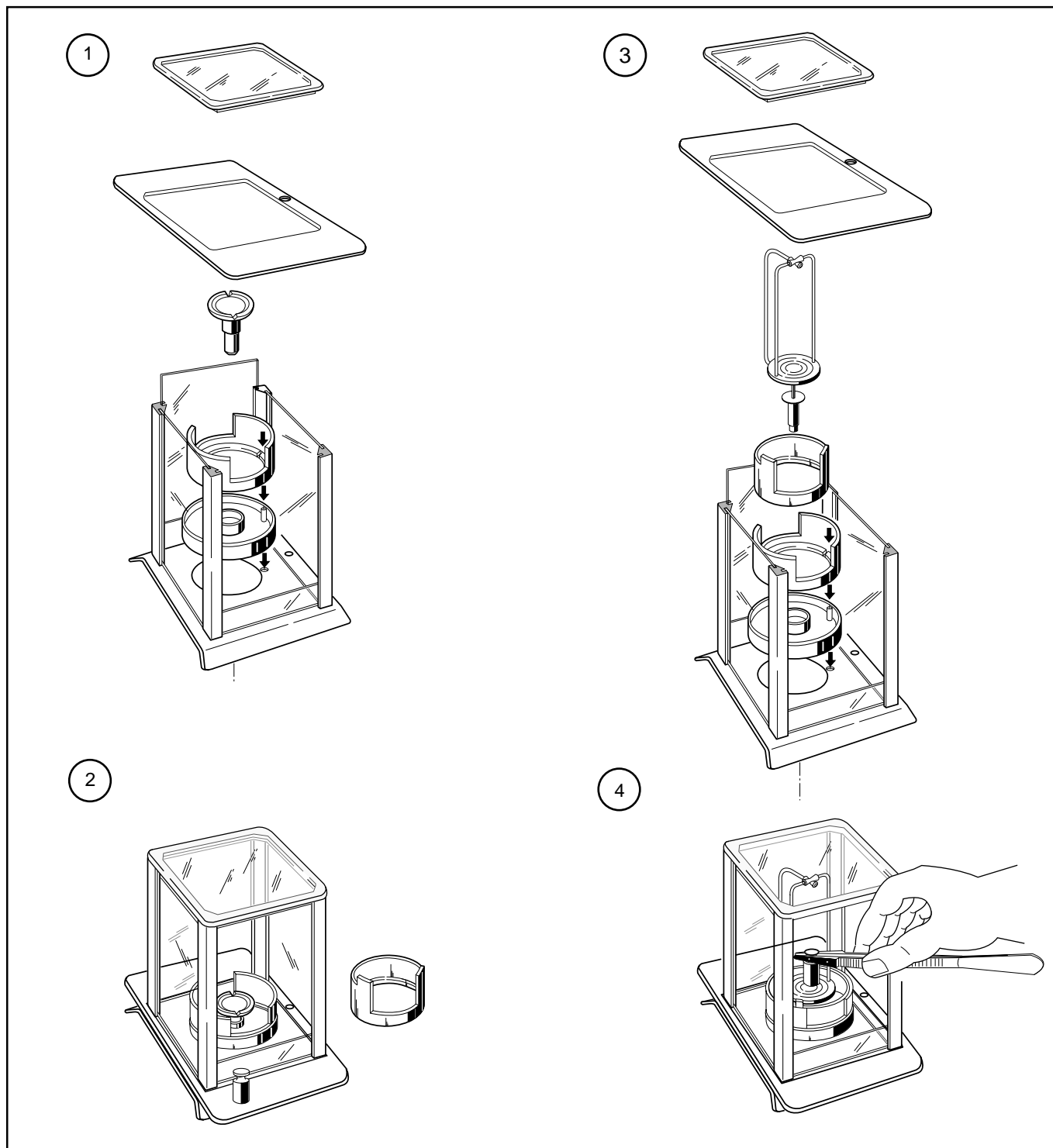

With the **METTLER TOLEDO AT21/AX26** with 1 microgram readability, the avoidance of shocks and air currents during the weighing process is of immense importance. This requirement is lent positive support by the evenly motorized door actuation; the motor coupling has thus been dispensed with. Possible manual door actuation is achievable by suitable configuration of the coupling elements, on the other hand the opening and closing of the right door with the left hand and vice versa is not possible here.

#### **Standard accessories:**

##### **Suspended balance pan**

The suspended pan prevents possible cornerload errors. If you nevertheless wish to work with the top loading weighing pan, you should ensure that the center of mass of the weights is in the middle of the weighing pan.

##### **Weight holder**

With the supplied weight holder, weights can be acclimatized in the weighing chamber. This prevents temperature differences between weighing sample and weighing chamber, which give rise to air currents.

Further tips regarding the proper way to work with electronic analytical and microbalances can be found in the enclosed **METTLER TOLEDO booklet "Weighing the Right Way"**.

##### **Default settings**

The balance has been configured for optimum operation under normal conditions. However, individual adjustments such as manual door operation, more restrictive stability detection (ASD), Auto-Zero off, etc. may be appropriate to achieve the best possible reproducibility.

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