RAININ

Pipetting 360°

Pipet Controller / Pipettierhelfer/
Accessoires de pipetage / Auxiliar de pipeteado

Pipet-X

Operating Manual
Gebrauchsanleitung
Mode d'emploi
Instrucciones de manejo

METTLER TOLEDO

Contents

Inhalt

1. Safety Instructions	4	1. Sicherheitsbestimmungen	4
2. Components	6	2. Einzelteile	6
3. Storage Conditions	7	3. Lagerbedingungen	7
4. Application	7	4. Funktion	7
5. Operating exclusions	7	5. Einsatzausschlüsse	7
6. Getting started	8	6. Die ersten Schritte	8
7. Pipetting settings	9	7. Pipettierfunktionen	9
8. Pipetting	10	8. Pipettieren	10
9. Recharging the battery	12	9. Akku laden	12
10. Replacing the battery	14	10. Akku austauschen	14
11. Battery disposal	16	11. Akkuentorgung	16
12. Cleaning	17	12. Reinigung	17
13. Replacing the filter and cleaning the check valve	17	13. Filter austauschen und Rückschlagventil reinigen	17
14. Autoclaving	19	14. Autoklavieren	19
15. Leak Test	20	15. Dichtigkeit prüfen	20
16. Troubleshooting	21	16. Störung - was tun?	21
17. Return for Repair	22	17. Zur Reparatur einsenden	22
18. Warranty	23	18. Mängelhaftung	23
19. Ordering Data	24	19. Bestelldaten	24
20. Accessories and spare parts	24	20. Zubehör und Ersatzteile	24

Sommaire

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1. Safety Instructions

Because this instrument may be used with hazardous materials, this Operating Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whomever uses this instrument to consult and establish appropriate safety and health practices and to determine the applicability of regulatory limitations prior to use.

1. Every user must read and understand this operating manual before operation.
2. Follow the general instructions for hazard prevention and safety regulations, e.g., wear protective clothing, eye protection and gloves.
3. Observe all specifications provided by reagent manufacturers.
4. Never use the instrument in an atmosphere with danger of explosion. Highly inflammable liquids must not be pipetted.
5. Use the instrument only for pipetting liquids, with strict regard to the defined limitations of use. If in doubt, contact the manufacturer or supplier.
6. Always use the instrument in such a way that neither the user nor any other person is endangered. Avoid splashes. Only dispense into suitable vessels.
7. Never use force on the instrument.
8. Use only original manufacturer's accessories and spare parts. Do not attempt to make any technical

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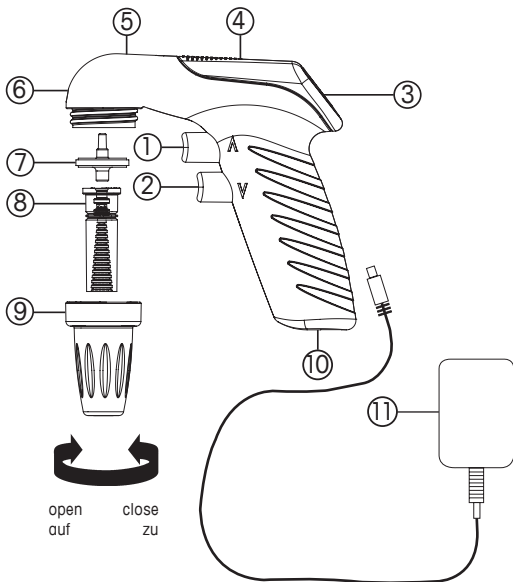
Warning!

Improper use of the instrument or the battery (short-circuit, mechanical damage, overheating, etc.) may cause the explosion of the battery.

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2. Components

1. Filling button
2. Dispensing button
3. Mode selection switch
4. Battery compartment
5. Recharging-indicator (LED)
6. Air inlet/exhaust
7. Membrane filter (PP/PTFE)
8. Adapter with check-valve (SI/PTFE)
9. Adapter support (PP)
10. Charging socket
11. AC adapter:
Input: AC 100 - 240 V; 50/60 Hz
Output: DC 5 V/200 mA

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3. Storage conditions

Store instrument and accessories in a cool, dry place.

Storage temperature:
- 20 °C to + 50 °C
(- 4 °F to + 122 °F)

4. Application

The instrument is designed to assist the filling and dispensing of graduated and volumetric pipettes of glass or plastic in the volume range of 0.1 ml to 100 ml for measuring liquids.

If the instrument is used correctly, the pipetted liquid will only contact the pipette. The instrument is not designed for Pasteur pipettes.

5. Operating exclusions

Never use the instrument with liquids whose vapors attack the materials silicone or EPDM.

Working- and recharging temperature:
+ 10 °C to + 35 °C
(+ 50 °F to + 95 °F)

Warning!

Never use or recharge the instrument in an atmosphere with danger of explosion. Highly inflammable liquids (e. g., ether, acetone and other liquids with a flash point below 0 °C) must not be pipetted.

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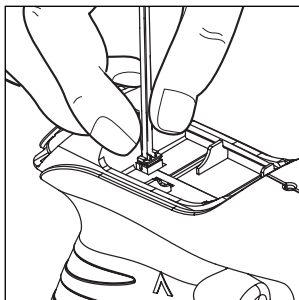
6. Getting started

1. Confirm alignment of battery plug and grip the plug by the wire and push it firmly into the socket. Place battery in compartment.

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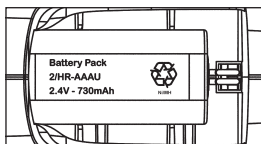
Note:

Flashing LED indicates that the battery has to be charged for initial operation (> pages 12-13).

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
2. Take battery compartment cover out of the packaging.

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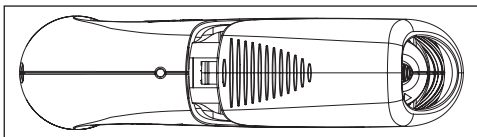▼ Blow-out
(with motor power)

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
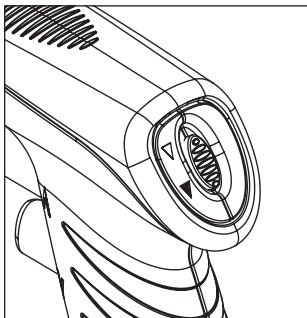
Pipetting speed control

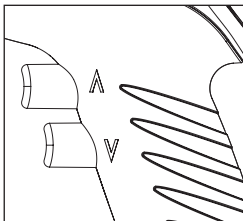The amount of pressure on the pipetting buttons provides continuously variable speeds up to the rate set by the speed controller. Press upper pipetting button for filling and lower button for dispensing.

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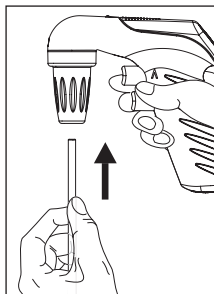
8. Pipetting

Attach the pipette

Hold the pipette as near to its upper end as possible, and carefully insert it into the adapter until it fits tightly.

Warning!


Be sure that the pipette fits tightly into the adapter. Never use force. Thin pipets are particularly liable to break. Avoid the risk of injury!

Once the pipette has been securely attached, always hold the instrument in a vertical position, tip down.

Filling

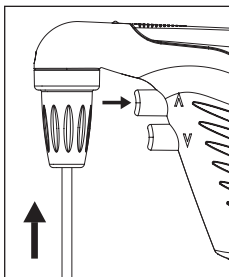Immerse the pipette tip into the liquid. Slowly depress the upper pipetting button and fill the pipette so that the meniscus is slightly above the mark desired.

Attention!

Take care to avoid overfilling of the pipette.

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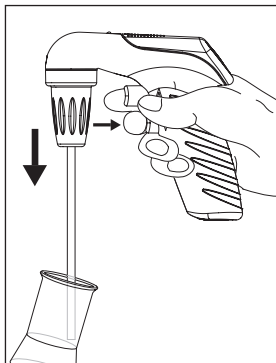
Adjust the volume

Use suitable lint-free tissue to wipe the pipette tip. Slowly press the lower button and dispense liquid until the meniscus is adjusted exactly to the desired volume.

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Dispensing

Slowly press the lower pipetting button. The dispensing speed is dependent on the amount of pressure on the pipetting button.

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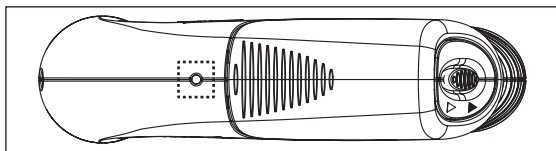
9. Recharging the battery

Warning!

Use original manufacturer's AC adapter only! Using a different AC adapter can damage both the instrument and the AC adapter, and will void the warranty. Confirm the voltage of the charger matches the voltage of your electrical outlet. The power supplied to the AC adapter should differ by no more than $\pm 10\%$.

One full charge of the nickel-metal hydride battery allows 8 hours of non-stop pipetting. A flashing LED indicates the need for recharging.

Recharging indicator

LED off or flashes green:

- battery charge is sufficient

LED flashes slowly red:

- remaining battery charge is sufficient for a maximum of two hours of pipetting; charging of battery is recommended

LED flashes quickly red:

- finish pipetting activity and charge battery

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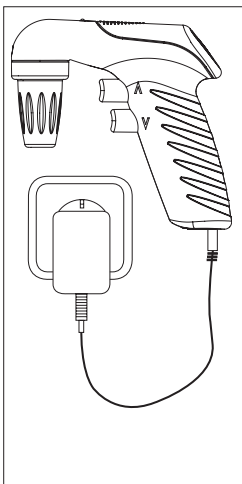
Recharging procedure

1. Temperature range for recharging:
+ 10 °C to + 35 °C
(+ 50 °F to + 95 °F)
2. When the AC adapter is initially connected, the LED on the instrument lights continuously.
3. After 4 hours of charging, the AC adapter switches to a pulsed, long-term charging mode (indicated by the LED flashing every 15 seconds).

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Note:

- The charging system is designed to prevent the battery from overcharging and minimizes the lazy-battery-effect. To maintain maximum battery capacity, it is best to charge the battery only when the slowly-flashing or quickly-flashing LED indicates charging is needed.
- The instrument can be used during recharging.
- If the LED does not light up during charging see chapter 'Troubleshooting' (➤ page 21).

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10. Replacing the battery

Warning!

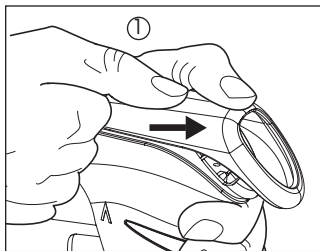
Only use original manufacturer's accessories (page 24). The battery must not be replaced by non-rechargeable or rechargeable batteries of other manufacturers. Use of the wrong batteries or improper use of the instrument (e.g., short-circuit, mechanical damage, overheating, etc.) may cause the batteries to explode.

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

- To open the battery compartment press at the embossed arrow, position ①, and simultaneously slide cover back.

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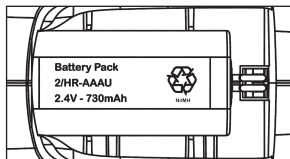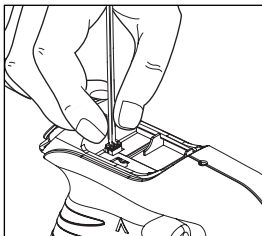11. Battery Disposal

The adjoining symbol means that storage batteries and electronic devices must be disposed of separately from household trash (mixed municipal waste) at the end of their service life.

- According to the Directive 2002/96/EC of the European Parliament and of the Council on Waste Electrical and Electronic Equipment (WEEE) of 27 January 2003, electronic equipment requires disposal according to the relevant national disposal regulations.
- Batteries contain substances that can have harmful effects on the environment and human health. Therefore according to the Directive 2006/66/EC of the European Parliament and the Council on Waste Batteries of 6 September 2006 batteries require disposal according to the relevant national disposal regulations. Dispose of batteries only when completely discharged.

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Warning!

Do not short-circuit the battery to discharge it!

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12. Cleaning

Occasionally wipe the housing with a damp cloth. When properly used, the instrument is maintenance-free.

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13. Replacing the filter and cleaning the check valve

Only necessary if a pipette has been overfilled or the suction capacity is impaired.

Warning!

Wear protective gloves and eye protection. Avoid the risk of contamination!

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
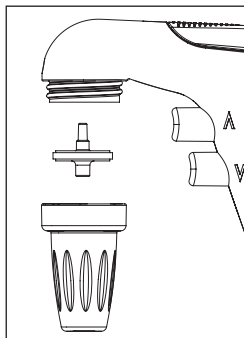
1. Pull out the pipette.
2. Unscrew the adapter support.
3. Pull out the membrane filter.

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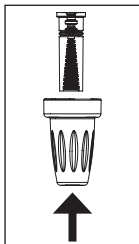
4. Remove the adapter from the adapter support by pushing it upwards. Do not remove check valve from the adapter.

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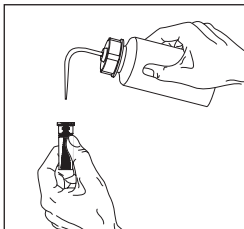5. To clean adapter with built-in check valve, carefully rinse it by using a wash bottle. Blow out the liquid and allow to dry completely.

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
6. Place the new membrane filter with the thicker end facing downwards into the adapter (with the smaller conical side up).

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14. Autoclaving

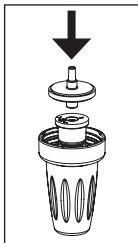1. The entire pipette adapter incl. membrane filter is autoclavable at 121°C (249,8 °F) (2 bar) according to DIN EN 285.

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Note:

Not every membrane filter is suitable. Only use genuine manufacturer's recommended accessories.

Each instrument will be supplied with a 0.2 µm membrane filter (red color code). The membrane filter is autoclavable up to 5 times at 121 °C (249,8 °F). It is the user's responsibility to ensure effective autoclaving. The filter with 0.2 µm pore size (for cell culture) is sufficient to meet class B tolerances. For class A accuracy, it is necessary to use a 3 µm membrane filter. This filter has a larger pore size and thus permits better airflow for gravity-delivery (Spare parts page 24).

2. Reassemble the instrument in reverse order and carry out leak test.

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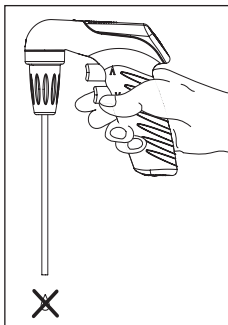
15. Leak Test

When the meniscus has been set, liquid should not drip out of the pipette until the valves are activated.

Should the pipette drip, see chapter 'Trouble-shooting' (> page 21)

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16. Troubleshooting / Störung – was tun?

Trouble	Possible Cause	Action to be taken
Suction capacity impaired	<ul style="list-style-type: none"> - Check valve closed - Filter or check valve dirty 	<ul style="list-style-type: none"> - Set mode selection switch for "power delivery" and press lower button for blow out (> p. 10) - Clean and/or replace check valve (> p. 17-20)
Pipette drips	<ul style="list-style-type: none"> - Filter not properly positioned - Adapter or pipette damaged 	<ul style="list-style-type: none"> - Insert filter properly (> p. 17) - Replace adapter or pipette (> p. 17-20)
Pipette not held properly	<ul style="list-style-type: none"> - Adapter dirty or damaged 	<ul style="list-style-type: none"> - Clean the adapter, allow it to dry and/or replace it (> p. 17-20)
After connecting the charger, the LED-display of the instrument flashes or does <u>not</u> light up	<ul style="list-style-type: none"> - Battery/Charger or Instrument defective 	<ul style="list-style-type: none"> - Replace battery/charger - Send the instrument in for repair (> p.22)
Recharging procedure not successful, LED-display of the instrument does not light up	<ul style="list-style-type: none"> - Battery problem 	<ul style="list-style-type: none"> - Replace battery (> p. 14, 15)

If recommendations in the Troubleshooting section do not solve current problems, contact your authorized dealer or the manufacturer directly.

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17. Return for Repair

Attention!

Transporting of hazardous materials without a permit is a violation of federal law.

RAININ will not accept instruments that are not appropriately cleaned and decontaminated.

Therefore, contact RAININ and obtain return authorization before sending your instrument for service.

Return the instrument, with the Return Authorization Number prominently displayed on the outside of the package to the address provided with the Return Authorization Number. Include an exact description of the type of malfunction and the media used.

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18. Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the instrument or the consequences of normal wear and tear especially of wearing parts such as pistons, seals, valves and the breakage of glass as well as the failure to follow the instructions of the operating manual. We are not liable for damage resulting from any actions not described in the operating manual or if non-original parts have been used.

Subject to technical modification without notice. Errors excepted.

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19. Ordering Data

Pipet-X - pipet controller with AC adapter,

incl. battery, battery compartment cover, 2 spare membrane filters (0.2 µm) and operating manual

with AC adapter for	Cat. No./MT Ordering
USA 120 V/60 Hz	PX-100/17011733
Europe (continental) 230 V/50 Hz	PX-100EU/17011743

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